

WARNING NOTICE: The experiments described in these materials are potentially hazardous and require a high level of safety training, special facilities and equipment, and supervision by appropriate individuals. You bear the sole responsibility, liability, and risk for the implementation of such safety procedures and measures. MIT shall have no responsibility, liability, or risk for the content or implementation of any of the material presented. [Legal Notices](#)

2. Purification of Solids by Recrystallization

2.1. Competent Chemist Rating: “How do you recrystallize a Mothball?”

Techniques Checklist:

- Solubility tests
- Choosing a good solvent system
- Inducing crystallization
- Filtration

Pre-Lab Discussion and Required Reading:

- Theory of recrystallization : Zubrick Ch. 13, LLP Ch. 11.2

Digital Lab Techniques Manual:


- 9. Recrystallization

Equipment:

- Test tubes (five 13x100 mm)
- Erlenmeyer flasks (2x50-mL, 1x125-mL)
- Small magnetic stir bar
- Funnel
- Filter paper
- Büchner funnel and filter paper
- Magnetic stirring/hot plate
- Filter flask (250-mL)
- Rubber filter adapter
- Large vial with white cap
- Test tube rack
- Large crystallizing dish
- Desiccator

Goal:

You will be given 2.00 g of impure naphthalene (mothballs!). Your job is to purify the naphthalene by recrystallization without losing a significant amount of your sample!¹

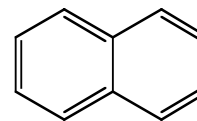

Naphthalene

Experiment Outline:

I. Solubility Tests

- Determine an appropriate solvent system for the recrystallization of naphthalene. For your tests try: **water, methanol, acetone, hexane, and toluene.** To understand how to find the appropriate solvent or solvent mixture for recrystallization, see pages 104–105 in Zubrick.

II. Recrystallization of Naphthalene

- Transfer the material to a 50-mL Erlenmeyer flask equipped with a stir bar. Add about 20 mL of the solvent (determined in Part I) and heat to boiling on a stir/hot plate.
- Remove any insoluble impurities by filtration, and recrystallize your product - *see Two-Solvent Recrystallization Guide.*
- Collect your crystals on a small Büchner funnel by vacuum filtration, and rinse with the cold solvent mixture.
- Your crystals should be colorless. If some orange or yellow color persists, wash your material with cold hexane. (*Be careful: What is the solubility of naphthalene in hexane?*).
- Dry your compound well - *see Two-Solvent Recrystallization Guide for tips.*
- Determine the yield, and obtain a melting point.

Results:

- To obtain your "CC Rating" in Purification of Solids by Crystallization, you must obtain colorless, well dried crystals weighing at least 1.30 g (no traces of yellow!). The purified material must melt over no more than three degrees with the lower range beginning no lower than 77 °C and the upper range ending no higher than 83 °C. This material must also be submitted to the TA for possible weight and melting point verification.

¹Adapted from Fieser, L. F.; Williamson, K. L. *Organic Experiments*; 7th ed.; D. C. Heath and Company: Lexington, MA, 1992; p. 40.